



## Molecular Devices Paradigm Plate Reader Configuration for Valita®Titer

Parameter	Instrument Settings	
Product	Valita®Titer (#VAL003)	Valita®Titer Plus (#VAL004)
Plate Type	96 Well Corning Half Area opaque	96 Well Corning Half Area opaque
Measurement Mode	Fluorescence Polarization	Fluorescence Polarization
Excitation	485 nm	485 nm
Emission	535 nm	535 nm
G Factor	1.0	1.0
Integration time	400 ms	400 ms
Read Height (mm)	4.1	4.8